

## Tomato + Minerals

Medium Detail Information Medium No.029MediumTomato + MineralsCompositionTomato 200g , grinded , boiled in 15-20 minutes and filtered before use

Glucose	10g/l
Yeast extract	10g/l
CaCO <sub>3</sub>	2g/l
Solution A	2.5ml/l
Solution B	2.5ml/l
Distilled water	1L

pH 6.9 ± 0.2 at 25°C

Solution A:	KH <sub>2</sub> PO <sub>4</sub>	10g/l
	K <sub>2</sub> HPO <sub>4</sub>	10g/l
	Distilled water	1 L
Solution B:	MgSO <sub>4</sub> .7 H <sub>2</sub> O	40g/l
	NaCL	2g/l
	FeSO <sub>4</sub>	2g/l
	MnSO <sub>4</sub>	2g/l
	Distilled water	1 L Comment